

Instructions Manual

ViewOne™ LabLite



ViewOne™ LabLite Instructions

1. Plug the LabLite into the receptacle on the universal AC adapter (Model: 3A-061WP05, output= 5V, 1.2A, input 100-240V, ~50-60Hz).
2. Plug the adapter into the power source.
3. Press the “Power” button once to turn on the LabLite and toggle to adjust intensity (high, medium, low)
4. Pressing the button while on low intensity will turn off the Lablite.



Note:

- The universal AC adapter comes with 4 different plug configurations. Before using, you must install the appropriate plug corresponding to the outlet of your country.
- Never attempt to open a ViewOne™ LabLite: This will compromise its function and will void any warranty.

I want another one!!!



Call, fax or email to tell us how many you want...

| Catalog # | Description | DIM |
|-----------|--------------------------|--------------------------------|
| VI - 1000 | ViewOne Lablite 100-240v | 17 (l) x 11.5 (w) x 0.6 (h) cm |

ViewOne Lablite

Order One Today...
View One Tomorrow!



7738 Arjons Drive • San Diego, CA 92126 • 800-255-1777 • 858 684-3190 • 858 684-3195 (fax) • inquiry@embitec.com
Embi Tec® is a registered trademark of C.C. IMEX. ViewOne™ is a trademark of Embi Tec

Related Products!

Use Tracking and Start Relaxing...



Pipetting a 96-well plate forcing you into solitary confinement? The LightOne™ Illuminator offers you the freedom to stop, work on something else and resume pipetting with full confidence of where you are. With the LightOne Illuminator, that 96-well plate should have ever been something you dreaded.



7738 Arjons Drive • San Diego, CA 92126 • 800-255-1777 • 858 684-3190 • 858 684-3195 (fax) • inquiry@embitec.com

Embi Tec® is a registered trademark of C.C. IMEX. ViewOne™ is a trademark of Embi Tec